

Mercury Collector, BÜRKLE

Bürkle's mercury collector enables the easy collection of spilt mercury, which can then be used again.

The mercury collector is composed of a PE tub. Within the screw cap lid there is a special plastic foam insert. On pressing this foam gently on the spilt mercury, the foam's pores are opened and the mercury globules are absorbed. When raised, the foam's pores are sealed again, thus keeping hold of the mercury.

The cap is then screwed on the tub, the globules are released again and fall through the perforated separation into the tub.



WWR4Q056



Article	Order No.
① Mercury collector	10 04 58901

