

## Pipette Controller VITLAB maneus®, VITLAB

The VITLAB maneus® pipette controller enables both left- and right-handers to operate all normal bulb and graduated pipettes from 0,1 to 200 ml easily and fatigue-free. Its safe and easy handling allows even inexperienced users to adjust the meniscus precisely.

With the new design, unscrewing the adapter enables easy and fast replacement of the hydrophobic membrane filter, which protects the instrument against fluid penetration.

The valve system is optimised so that liquids can be drawn up simply, without exerting pressure. The highly sensitive filling and discharge of liquids are controlled gently by the pipetting knob. Thus, the suction element provides rapid filling of the pipette (capacity: 50 ml in less than 10 seconds). The discharge bellows are used for the emptying (blow-out) of the pipette. The specially moulded intake cone ensures secure seating for all normal bulb and graduated pipettes (0.1 to 200 ml).

The VITLAB maneus® is simple to dismantle, easy to clean, and completely autoclavable at 121°C (2 bar) according to DIN EN 285.

With replacement 3 µm membrane filter and instruction manual.

| Article                           | Order No.   |
|-----------------------------------|-------------|
| Pipette controller VITLAB maneus® | 10 85 30500 |

